

**IMPORTANT WARNING WHEN REPLACING A CONTROLLER ON OPxxx, OMAXI, OMINI**

On some old osmosers, the cable bundle has no white wire on the 4th pin of the connector.



old cable bundle:

standard OPxxx, OMAXI, OMINI cable bundle:



---

**ANY CABLE BUNDLE WITHOUT WHITE WIRE ON PIN 4 MUST BE REPLACED BEFORE CONNECTING A NEW CONTROLLER !!!**

( in such a case, the internal fuse will burn and the osmoser will not work )

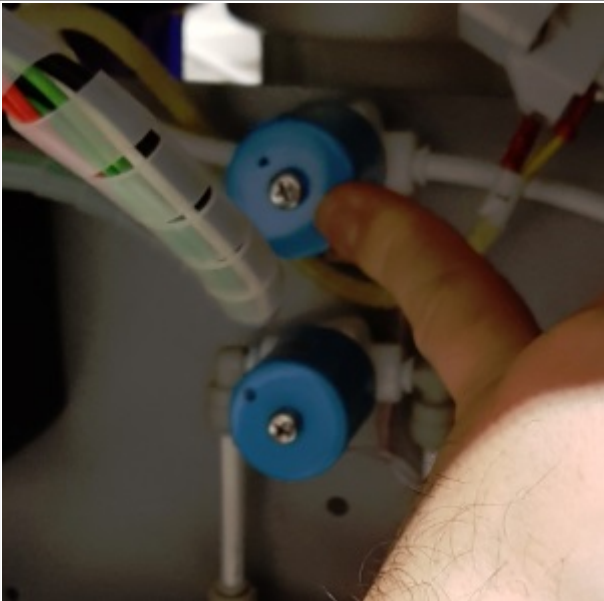
---

## CONNECTION OF THE ELECTRIC CABLE BUNDLE

[Original file](#)



| WIRE                  | CONNECTED TO                                                                                                                        |
|-----------------------|-------------------------------------------------------------------------------------------------------------------------------------|
| <b>Cable 1 Yellow</b> | <br>Low Pressure Switch                          |
| <b>Cable 2 Blue</b>   | <br>Cable connector of the High Pressure Switch |

|                            |                                                                                                                |
|----------------------------|----------------------------------------------------------------------------------------------------------------|
| <b>Cable 3 Red</b>         |  <p>Low Solenoid Valve</p>    |
| <b>Cable 4 Green Black</b> | <div><div><div>✖</div><div>✖</div></div><p>T jonction</p></div>                                                |
| <b>Cable 5 Green White</b> |  <p>High Solenoid Valve</p> |

**Cable 6 Blue Brown**



24V Power supply

From:

<https://diasys-technologies.com/waterpurifiers/> - **DiaSys Water Purifier Systems**

Permanent link:

[https://diasys-technologies.com/waterpurifiers/doku.php?id=replace\\_controllerwithoutwhitewire&rev=1573833428](https://diasys-technologies.com/waterpurifiers/doku.php?id=replace_controllerwithoutwhitewire&rev=1573833428)



Last update: **2019/11/15 15:57**