

- verify the main power is active - RED Light in the power switch button must be ON. - verify the 24V output of the adaptor. Replace the adaptor if necessary.

- if 24VDC is OK; verify the fuse inside the controller box. **Before opening, switch off first the system.**

From:

<https://diasys-technologies.com/waterpurifiers/> - **DiaSys Water Purifier Systems**

Permanent link:

<https://diasys-technologies.com/waterpurifiers/doku.php?id=allindicatoroff&rev=1552903257> 

Last update: **2019/03/18 10:00**