


IMPORTANT WARNING WHEN REPLACING A CONTROLLER ON OPxxx, OMAXI, OMINI

On some old osmosers, the cable bundle has no white wire on the 4th pin of the connector.



old cable bundle:

standard OPxxx, OMAXI, OMINI cable bundle:

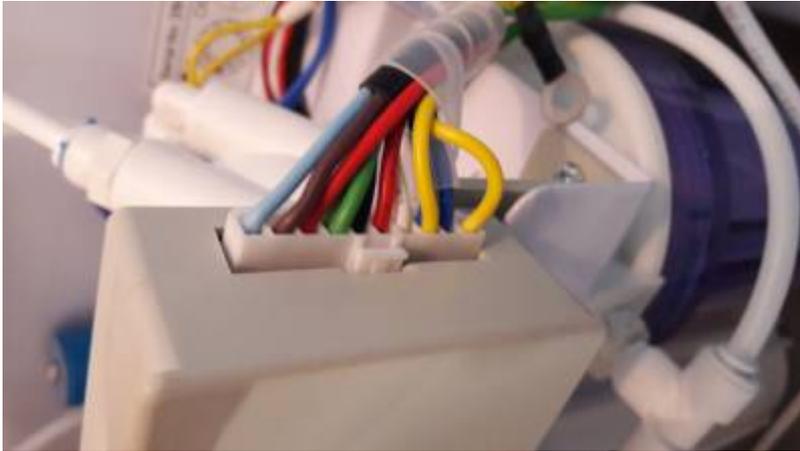

ANY CABLE BUNDLE WITHOUT WHITE WIRE ON PIN 4 MUST BE REPLACED BEFORE CONNECTING A NEW CONTROLLER !!! (in such a case, the internal fuse will burn and the osmoser will not work)

From:

<https://diasys-technologies.com/waterpurifiers/> - **DiaSys Water Purifier Systems**

Permanent link:

https://diasys-technologies.com/waterpurifiers/doku.php?id=replace_controllerwithoutwhitewire&rev=1573724221

Last update: **2019/11/14 09:37**