

- verify the main power is active - RED Light on the power switch button must be ON.
- verify the 24V output of the adaptor. Replace the adaptor if necessary.
- if 24VDC is OK; verify the fuse inside the controller box. **Before opening, first of all, switch off the system!**

From:

<http://diasys-technologies.com/waterpurifiers/> - **DiaSys Water Purifier Systems**

Permanent link:

<http://diasys-technologies.com/waterpurifiers/doku.php?id=allindicatoroff>

Last update: **2019/03/25 17:05**

