

to disconnect the internal autofilling system implemented in the OMINI-Autofilling version, proceed as follow:

1. unplug AC power supply
2. open the osmoser
3. remove the tubing connected to the purified water output
4. remove the tubing at the left of the bottom blue valve
5. connect the purified water output directly to the left of the bottom blue valve
6. close the osmoser
7. connect a hand valve on the purified water output
8. disconnect the float sensor cable
9. reconnect AC power supply
10. switch on the osmoser

From:

<http://diasys-technologies.com/waterpurifiers/> - **DiaSys Water Purifier Systems**

Permanent link:

<http://diasys-technologies.com/waterpurifiers/doku.php?id=disconnect-autofilling>

Last update: **2019/03/15 15:34**

